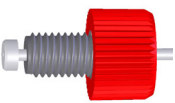
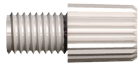
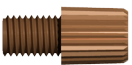
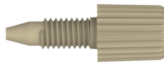



No Guesswork & Reusable!

Click-N-Seal® offers a unique system for hand-tightened fittings for flared or ferrule connections with the unique torque-control system. As the fitting is tightened, the resistance of the cap increases. When the correct torque is reached the cap clicks. The user is able to hand-tighten the fittings without tools or guesswork and be confident that the fitting is properly seated, creating a secure seal. The fitting can be unscrewed and reused.

- Available in Polycarbonate, Acetal or PEEK
- Standard or compact head for space saving design
- Compatible with Dibafit™ Type P and Type S ferrules or inverted cones
- Pressure rated up to 1000 psi (Click-N-Seal® Ultra 10-32 pressure rated up to 2000psi)
- Retains 1/16", 1/8" or 3/16" OD semi-rigid tubing e.g. PTFE, ETFE, FEP
- For 1/4"-28 UNF, M6, 10-32 or 6-40 ports

Click-N-Seal® Line *(shown at actual size)*

|  |  |  |  |  |
|---|---|--|--|---|
| Click-N-Seal® Regular | Click-N-Seal® Plus | Click-N-Seal® Ultra | Click-N-Seal® Ultra10-32 Cone Tip | Click-N-Seal® Micro 6-40 |
| <ul style="list-style-type: none"> • Eliminates over-and under-torquing • No tools required • Made of polycarbonate & acetal • Suitable for flat-bottom 1/4"-28 UNF or M6 ports | <ul style="list-style-type: none"> • Eliminates over-and under-torquing • No tools required • Made of acetal • Slim profile • Suitable for flat-bottom 1/4"-28 UNF or M6 ports | <ul style="list-style-type: none"> • Eliminates over-and under-torquing • No tools required • Made of PEEK • Slim profile • Works well with aggressive chemistries • Suitable for flat-bottom 1/4"-28 UNF, M6 or 10-32 ports | <ul style="list-style-type: none"> • Eliminates over-and under-torquing • No tools required • Made of PEEK • Slim profile • Works well with aggressive chemistries • Suitable for HPLC type ports • Also available in flat bottom | <ul style="list-style-type: none"> • Eliminates over-and under-torquing • No tools required • Made of PEEK • Slim profile • Works well with aggressive chemistries • Suitable for flat-bottom 6-40 ports • Ideal for microfluidic applications |


Click-N-Seal® Micro 6-40 Fitting

The new Click-N-Seal® Micro 6-40 fitting provides torque control in the smallest spaces

NEW

Ordering Information

| CLICK-N-SEAL® MICRO FOR 1/16" OD TUBING / 6-40 THREAD | | | | |
|---|----------|---------|--------|-------|
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF16-KT9L | PEEK™ | Natural | 6-40 | 10 pk |
| 008NF16-KT9B | PEEK™ | Black | 6-40 | 10 pk |
| 008NF16-KT9R | PEEK™ | Red | 6-40 | 10 pk |

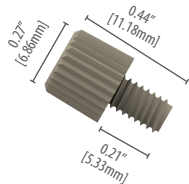


008NF16-KT9X
Click-N-Seal® Micro
6-40 fitting

10-32 Click-N-Seal® Ultra Fittings

The 10-32 Cone-Tip Click-N-Seal® fitting solution for quick and easy connection of 1/16" tube into a 10-32 conical port

Ordering information

| CLICK-N-SEAL® FITTINGS FOR 1/16" OD TUBING / 10-32 THREAD | | | | |
|---|----------|---------|--------|-------|
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NT16-KT4L | PEEK™ | Natural | 10-32 | 10 pk |

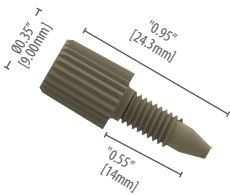


008NT16-KT4L
10-32 cone-tip Click-N-Seal®
fitting

The 10-32 flat bottom Click-N-Seal® fitting solution for quick and easy connection of 1/16" tube into a 10-32 flat bottom port

Ordering information

| CLICK-N-SEAL® FITTINGS FOR 1/16" OD TUBING / 10-32 THREAD | | | | |
|---|----------|---------|--------|-------|
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF16-KT4L | PEEK™ | Natural | 10-32 | 10 pk |

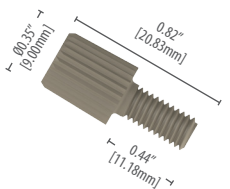


008NF16-KT4L
10-32 flat bottom Click-N-Seal®
fitting

Click-N-Seal® Fittings For Flares


Click-N-Seal® fittings for flared tubing enable you to create a ferrule-free connection and are offered in three different bore sizes - simply choose the size that best matches your tube OD. We recommend the use of a washer to ensure the flare seals correctly. An o-ring may also be necessary to ensure the fitting seats correctly in deeper ports and in this case a second washer should be used between the o-ring and the fitting.

Ordering information

| CLICK-N-SEAL® FITTINGS FOR FLARED TUBING UP TO 1/16" OD | | | | |
|--|----------|---------|--|-------|
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG16-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NG16-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NG16-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NG16-CS6B | PC | Black | M6 | 10 pk |
| 008NG16-CS6H | PC | White | M6 | 10 pk |
| 008NG16-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR FLARED TUBING UP TO 1/16" OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG16-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NG16-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR FLARED TUBING UP TO 1/16" OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG16-KT5L | PEEK™ | Natural | 1/4-28 | 10 pk |
| 008NG16-KT4L | PEEK™ | Natural | 10-32 | 10 pk |
| CLICK-N-SEAL® FITTINGS FOR FLARED TUBING 1.6MM TO 2.5MM OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG25-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NG25-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NG25-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NG25-CS6B | PC | Black | M6 | 10 pk |
| 008NG25-CS6H | PC | White | M6 | 10 pk |
| 008NG25-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR FLARED TUBING 1.6MM TO 2.5MM OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG25-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NG25-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR FLARED TUBING 1.6MM TO 2.5MM OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG25-KT5L | PEEK™ | Natural | 1/4-28 | 10 pk |
| CLICK-N-SEAL® FITTINGS FOR FLARED TUBING 2.5MM TO 1/8" OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG32-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NG32-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NG32-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NG32-CS6B | PC | Black | M6 | 10 pk |
| 008NG32-CS6H | PC | White | M6 | 10 pk |
| 008NG32-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR FLARED TUBING 2.5MM TO 1/8" OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG32-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NG32-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR FLARED TUBING 2.5MM TO 1/8" OD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NG32-KT5L | PEEK™ | Natural | 1/4-28 | 10 pk |
| WASHER & O-RING KITS | | | PART NUMBER | QTY |
| DESCRIPTION | | | | |
| 008FLK-16M | | | 2 x PEEK™ washer & 1 x FKMFFM O-ring for 1/16" | 10 pk |
| 008FLK-25M | | | 2 x PEEK™ washer & 1 x FKMFFM O-ring for 2.5MM | 10 pk |
| 008FLK-32M | | | 2 x PEEK™ washer & 1 x FKMFFM O-ring for 1/8" | 10 pk |

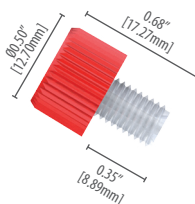

008NG16-CS/xx¹
Click-N-Seal® standard fitting
for 1/16" OD tubing

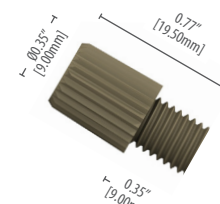

008NG16-ATxx¹/KT5L
Click-N-Seal® Plus/ Ultra fitting
for 1/16" OD tubing

¹The dimensions shown apply to all variants within the product group.

Click-N-Seal® Self-Torquing Adaptors

Thread to barb connections that can't be overtightened!


Ordering information

CLICK-N-SEAL® MALE 1/4"-28 UNF TO BARB ADAPTOR


| PART NUMBER | MATERIAL | COLOR | THREAD | TUBE ID | QTY |
|--------------|----------|-------|----------|----------------|-------|
| 008NB16-CS5B | PC | Black | 1/4•28 M | 1/16" (1.57mm) | 10 pk |
| 008NB16-CS5H | PC | White | 1/4•28 M | 1/16" (1.57mm) | 10 pk |
| 008NB16-CS5R | PC | Red | 1/4•28 M | 1/16" (1.57mm) | 10 pk |
| 008NB16-CS6B | PC | Black | M6 | 1/16" (1.57mm) | 10 pk |
| 008NB16-CS6H | PC | White | M6 | 1/16" (1.57mm) | 10 pk |
| 008NB16-CS6R | PC | Red | M6 | 1/16" (1.57mm) | 10 pk |
| 008NB32-CS5B | PC | Black | 1/4•28 M | 1/8" (3.2mm) | 10 pk |
| 008NB32-CS5H | PC | White | 1/4•28 M | 1/8" (3.2mm) | 10 pk |
| 008NB32-CS5R | PC | Red | 1/4•28 M | 1/8" (3.2mm) | 10 pk |
| 008NB32-CS6B | PC | Black | M6 | 1/8" (3.2mm) | 10 pk |
| 008NB32-CS6H | PC | White | M6 | 1/8" (3.2mm) | 10 pk |
| 008NB32-CS6R | PC | Red | M6 | 1/8" (3.2mm) | 10 pk |
| 008NB40-CS5B | PC | Black | 1/4•28 M | 3/32" (4.0mm) | 10 pk |
| 008NB40-CS5H | PC | White | 1/4•28 M | 3/32" (4.0mm) | 10 pk |
| 008NB40-CS5R | PC | Red | 1/4•28 M | 3/32" (4.0mm) | 10 pk |

CLICK-N-SEAL® ULTRA MALE 1/4"-28 UNF TO BARB ADAPTOR

| PART NUMBER | MATERIAL | COLOR | THREAD | TUBE ID | QTY |
|--------------|----------|---------|----------|----------------|-------|
| 008NB16-KT5L | PEEK™ | Natural | 1/4•28 M | 1/16" (1.57mm) | 10 pk |
| 008NB16-KT6L | PEEK™ | Natural | M6 | 1/16" (1.57mm) | 10 pk |
| 008NB32-KT5L | PEEK™ | Natural | 1/4•28 M | 1/8" (3.2mm) | 10 pk |
| 008NB32-KT6L | PEEK™ | Natural | M6 | 1/8" (3.2mm) | 10 pk |

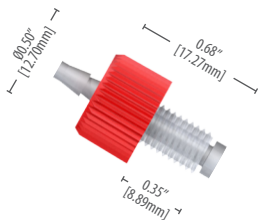

008NBxx¹-CS Series
Click-N-Seal® Thread to Barb

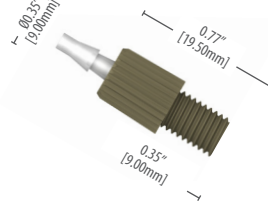

008NBxx¹-KT Series
Click-N-Seal® Ultra Thread to Barb

¹The dimensions shown apply to all variants within the product group. Barb diameters and lengths vary according to the ID of tube.

Dibafit™ Type P Ferrule System

Chemically inert flangeless fitting with permanently attached PTFE ferrule ideal for repeat connect/disconnect

Ordering information

FERRULES FOR 1/16" OD TUBING

| PART NUMBER | DESCRIPTION | QTY |
|-------------|---|-------|
| 008FT16 | PTFE / SS ferrule for 1/4•28 & M6 fittings | 10 pk |
| 008FTK16 | PTFE / PEEK™ ferrule for 1/4•28 & M6 fittings | 10 pk |

FERRULES FOR 1/8" OD TUBING

| PART NUMBER | DESCRIPTION | QTY |
|-------------|---|-------|
| 008FT32 | PTFE / SS ferrule for 1/4•28 & M6 fittings | 10 pk |
| 008FTK32 | PTFE / PEEK™ ferrule for 1/4•28 & M6 fittings | 10 pk |

CLICK-N-SEAL® FITTINGS FOR 1/16" OD TUBING

| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|-------|--------|-------|
| 008NF16-CS5B | PC | Black | 1/4•28 | 10 pk |
| 008NF16-CS5H | PC | White | 1/4•28 | 10 pk |
| 008NF16-CS5R | PC | Red | 1/4•28 | 10 pk |
| 008NF16-CS6B | PC | Black | M6 | 10 pk |
| 008NF16-CS6H | PC | White | M6 | 10 pk |
| 008NF16-CS6R | PC | Red | M6 | 10 pk |

CLICK-N-SEAL® FITTINGS FOR 1/8" OD TUBING

| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|-------|--------|-------|
| 008NF32-CS5B | PC | Black | 1/4•28 | 10 pk |
| 008NF32-CS5H | PC | White | 1/4•28 | 10 pk |
| 008NF32-CS5R | PC | Red | 1/4•28 | 10 pk |
| 008NF32-CS6B | PC | Black | M6 | 10 pk |
| 008NF32-CS6H | PC | White | M6 | 10 pk |
| 008NF32-CS6R | PC | Red | M6 | 10 pk |

CLICK-N-SEAL® PLUS FITTINGS FOR 1/16" OD TUBING

| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|-------|--------|-------|
| 008NF16-AT5H | Acetal | White | 1/4•28 | 10 pk |
| 008NF16-AT6H | Acetal | White | M6 | 10 pk |

CLICK-N-SEAL® PLUS FITTINGS FOR 1/8" OD TUBING

| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|-------|--------|-------|
| 008NF32-AT5H | Acetal | White | 1/4•28 | 10 pk |
| 008NF32-AT6H | Acetal | White | M6 | 10 pk |

CLICK-N-SEAL® ULTRA FITTINGS FOR 1/16" OD TUBING


| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|---------|--------|-------|
| 008NF16-KT5L | PEEK™ | Natural | 1/4•28 | 10 pk |

CLICK-N-SEAL® ULTRA FITTINGS FOR 1/8" OD TUBING


| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|---------|--------|-------|
| 008NF32-KT5L | PEEK™ | Natural | 1/4•28 | 10 pk |

BLANKING PLUG FOR 1/4•28 FLAT-BOTTOM PORTS

| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|-------------|----------|---------|--------|------|
| 002320 | ETFE | Natural | 1/4•28 | 2 pk |

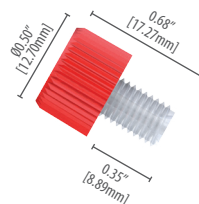

008NF16-CS/ATxx¹
Click-N-Seal® fitting
for 1/16" OD tubing

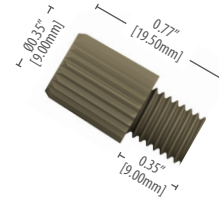

008NF16-ATxx¹/KT5L
Click-N-Seal® Plus/ Ultra fitting
for 1/16" OD tubing

¹The dimensions shown apply to all variants within the product group

Dibafit™ Gripper Fittings

Chemically inert flangeless fitting with permanently attached PTFE ferrule ideal for repeat connect/disconnect


Ordering information

GRIPPERS FOR 1/16" OD TUBING

| PART NUMBER | DESCRIPTION | QTY |
|-------------|---|-------|
| 002310 | PTFE / SS ferrule for 1/4•28 fittings | 10 pk |
| 002313 | PTFE / SS ferrule for M6 fitting fittings | 10 pk |
| 008FTK16 | PTFE / PEEK™ ferrule for 1/4•28 & M6 fittings | 10 pk |

CLICK-N-SEAL® ULTRA FITTINGS FOR 1/16" OD TUBING

| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
|--------------|----------|---------|--------|-------|
| 008NG16-KT5L | PEEK™ | Natural | 1/4•28 | 10 pk |
| 008NG16-KT4L | PEEK™ | Natural | 10-32 | 10 pk |

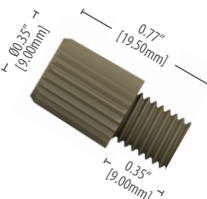

008NG16-KTXL
Click-N-Seal® Plus/ Ultra fitting
for 1/16" OD tubing

Dibafit™ Type S Ferrule System

Chemically inert with permanently attached PEEK™ or ETFE ferrules, ideal for repeated connect/disconnect

Dibafit™ Inverted Cone System

Removable and reusable system for quick and convenient low-pressure connections

Ordering information


| FERRULES FOR 1/16" OD TUBING | | | | |
|---|---|---------|--------|-------|
| PART NUMBER | DESCRIPTION | | | QTY |
| 008FK16 | PEEK™ ferrule and SS lock ring | | | 10 pk |
| 008FZ16 | ETFE ferrule and SS lock ring | | | 10 pk |
| 008FK16-C1 | PEEK™ ferrule; taper port and SS lock ring | | | 10 pk |
| FERRULES FOR 1/16" OD TUBING / 6-40 THREAD NEW | | | | |
| PART NUMBER | DESCRIPTION | | | QTY |
| 008FK16-03C9L | PEEK™ ferrule; taper port, 0.3mm through hole | | | 10 pk |
| 008FK16-05C9L | PEEK™ ferrule; taper port, 0.5mm through hole | | | 10 pk |
| 008FK16-08C9L | PEEK™ ferrule; taper port, 0.8mm through hole | | | 10 pk |
| 008FK16-F9L | PEEK™ ferrule; flat port, and SS lock ring | | | 10 pk |
| CLICK-N-SEAL® FITTINGS FOR 1/16" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF16-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NF16-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NF16-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NF16-CS6B | PC | Black | M6 | 10 pk |
| 008NF16-CS6H | PC | White | M6 | 10 pk |
| 008NF16-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR 1/16" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF16-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NF16-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR 1/16" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF16-KT4L | PEEK™ | Natural | 10-32 | 10 pk |
| 008NF16-KT5L | PEEK™ | Natural | 1/4-28 | 10 pk |
| CLICK-N-SEAL® MICRO FOR 1/16" OD TUBING / 6-40 THREAD | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF16-KT9L | PEEK™ | Natural | 6-40 | 10 pk |
| 008NF16-KT9B | PEEK™ | Black | 6-40 | 10 pk |
| 008NF16-KT9R | PEEK™ | Red | 6-40 | 10 pk |
| FERRULES FOR 1/8" OD TUBING | | | | |
| PART NUMBER | DESCRIPTION | | | QTY |
| 008FK32 | PEEK™ ferrule and SS lock ring | | | 10 pk |
| 008FZ32 | ETFE ferrule and SS lock ring | | | 10 pk |
| CLICK-N-SEAL® FITTINGS FOR 1/8" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF32-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NF32-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NF32-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NF32-CS6B | PC | Black | M6 | 10 pk |
| 008NF32-CS6H | PC | White | M6 | 10 pk |
| 008NF32-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR 1/8" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF32-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NF32-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR 1/8" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NF32-KT5L | PEEK™ | Natural | 1/4-28 | 10 pk |

Ordering information

| INVERTED CONES FOR 1/16" OD TUBING | | | | |
|---|--------------------|--------------|--------|-------|
| PART NUMBER | DESCRIPTION | | | QTY |
| 008CZ16 | ETFE inverted cone | | | 10 pk |
| CLICK-N-SEAL® FITTINGS FOR 1/16" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NC16-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NC16-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NC16-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NC16-CS6B | PC | Black | M6 | 10 pk |
| 008NC16-CS6H | PC | White | M6 | 10 pk |
| 008NC16-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR 1/16" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NC16-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NC16-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR 1/16" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NC16-KT5L | PEEK™ | Natural | 1/4-28 | 10 pk |
| INVERTED CONES FOR 2.0, 2.5, 3.0 & 3.2MM (1/8") OD TUBING | | | | |
| PART NUMBER | MATERIAL | TUBE SIZE | | QTY |
| 008CZ19 | ETFE | 1.8 - 2.0mm | | 10 pk |
| 008CZ25 | ETFE | 2.5mm | | 10 pk |
| 008CZ30 | ETFE | 3.0mm | | 10 pk |
| 008CZ32 | ETFE | 1/8" (3.2mm) | | 10 pk |
| 001545 | ETFE | 1/8" (3.2mm) | | 10 pk |

All of the cones listed above are compatible with any of the fittings also listed.


| CLICK-N-SEAL® FITTINGS FOR 1/8" OD TUBING | | | | |
|---|----------|---------|--------|-------|
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NC32-CS5B | PC | Black | 1/4-28 | 10 pk |
| 008NC32-CS5H | PC | White | 1/4-28 | 10 pk |
| 008NC32-CS5R | PC | Red | 1/4-28 | 10 pk |
| 008NC32-CS6B | PC | Black | M6 | 10 pk |
| 008NC32-CS6H | PC | White | M6 | 10 pk |
| 008NC32-CS6R | PC | Red | M6 | 10 pk |
| CLICK-N-SEAL® PLUS FITTINGS FOR 1/8" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NC32-AT5H | Acetal | White | 1/4-28 | 10 pk |
| 008NC32-AT6H | Acetal | White | M6 | 10 pk |
| CLICK-N-SEAL® ULTRA FITTINGS FOR 1/8" OD TUBING | | | | |
| PART NUMBER | MATERIAL | COLOR | THREAD | QTY |
| 008NC32-KT5L | PEEK | Natural | 1/4-28 | 10 pk |

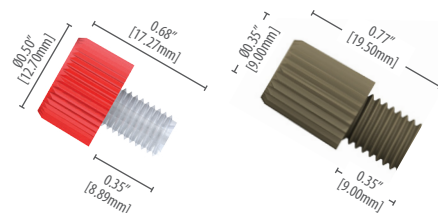

008NF16-CSxx¹
Click-N-Seal fitting for 1/16" OD tubing

008NF16-ATxx¹/KT5L
Click-N-Seal Plus/ Ultra fitting for 1/16" OD tubing

¹The dimensions shown apply to all variants within the product group.

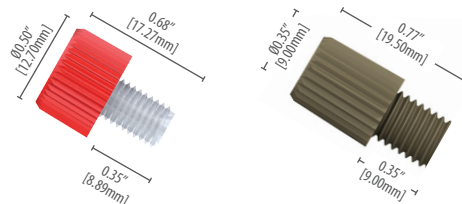

008NC16-CSxx¹
Click-N-Seal Fitting for 1/8" OD tubing

008NC16-ATxx¹/KT5L
Click-N-Seal Plus/ Ultra Fitting for 1/8" OD tubing

¹The dimensions shown apply to all variants within the product group.



United States
Diba Industries Inc.
4 Precision Road
Danbury, CT 06810 USA
Tel: +1 (203)-744-0773
Fax: +1 (203)-744-0663
salesdept@dibaind.com

Europe
Diba Industries Ltd.
2 College Park
Coldhams Lane
Cambridge, CB1 3HD, UK
Tel: +44 (0) 1223 472801
Fax: +44 (0) 1223 416787
sales@dibaind.com

China
Diba Industries Inc.
Halma International Ltd.
Shanghai Representative Office
Room 1106, Metro Plaza, No555
Lou Shan Guan Road,
Shanghai, PRC 200051, CHINA
Tel: +86-21-6016-7622
Fax: +86-21-5206-8191
chinasales@dibaind.com